

Quick Start Guide



Trigno® Avanti Mobile


NEED ASSISTANCE?


	SUPPORT	TRAINING
North America	+1 (508) 545-8200 support@delsys.com	+1 (508) 545-8200 training@delsys.com
Europe	+44 161 504 5066 support@delsyseurope.com	+44 161 504 5066 training@delsyseurope.com




1 Experiment Set-up

Follow steps **A** to **D** to achieve high quality surface EMG signals.


- A** Remove sensors from charging cradle and touch 

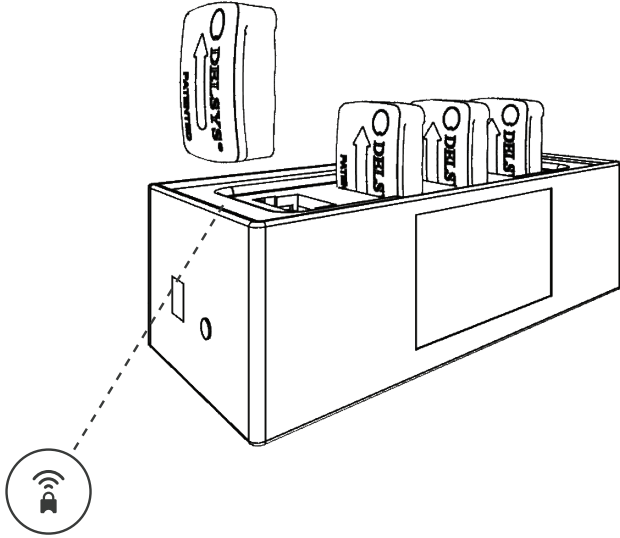


- B** Open the desired application from the Mobile EMG Suite in your Android tablet. App will automatically start scanning for sensors.

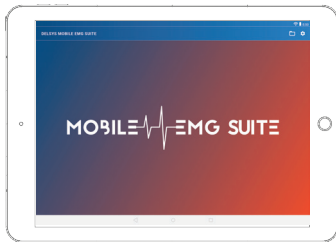

- C** Clean the shaved muscle sites with an alcohol wipe.

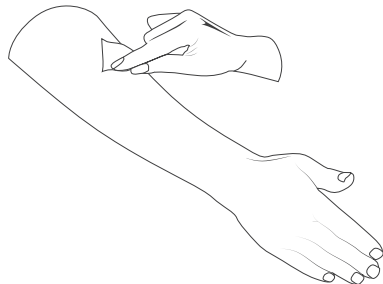




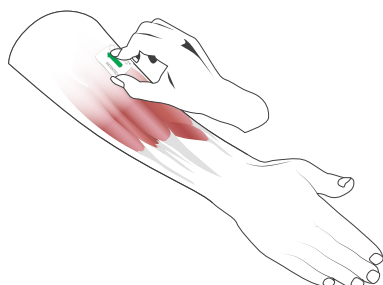

- D** Place the sensor(s) over the middle of the muscle(s) using Adhesive Interfaces.




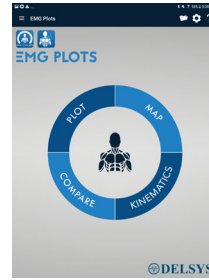

2 Record Data with Mobile EMG Suite

- A** Select your preferred acquisition module.



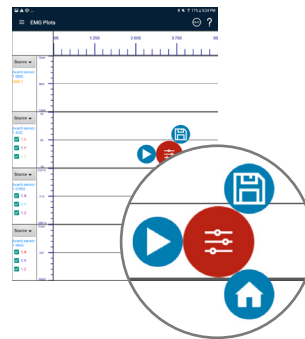


- B** Click  to show Start , Record , and Home  buttons.





- C** Click  to start data stream. Click  to record data.

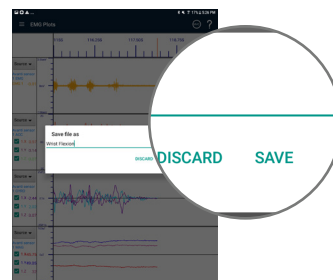

- D** Click  to stop recording. Edit the file name as desired and click **SAVE** to save data. Click Stop  to end data stream.



- E** Click the  icon in the top right, find your file, click , and select your preferred option for data management and export.

